Microscope Eyepiece Digital Cameras

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levenhuk 
Zoom&Joy®

Description

Levenhuk C NG Series eyepiece digital cameras have been designed exclusively for microscopes. They can be installed on any type of optical microscopes, be it biological or metallurgical, mono or stereo microscope. The observed view may be transferred directly to your PC, and with the software included in the kit you can view and edit the resulting image.

Features

- Eyepiece digital cameras for microscopes may be used with any type of optical microscopes. Whether you have a biological or a metallurgical microscope, whether it is mono or stereo, you can install the camera and use it straight away.
- C NG series includes cameras with resolutions ranging from 0.3 to 14.0 megapixels.
- The image can be transferred and saved on your PC exactly the way you see it in the oculars.
- The standard kit includes Levenhuk ToupView software necessary to view and edit your images which supports *.bmp, *.jpg, *.jpeg, *.png, *.tif, *.tiff, *.gif, *.psd, *.ico, *.emf, *.wmf, and other formats.
- The camera can be charged and connected to your PC with USB-cable.
- Compatible OS: Mac OS, Linux, Windows XP/Vista/7/8 (32-bit or 64-bit).

Package

- Digital camera
- USB cable
- 2 adapters for different microscope types
- Installation CD (setup wizard, drivers, Levenhuk ToupView software for image viewing and editing)
- User manual and warranty

Caution

- **Never look at the sources of bright light or lasers through the camera or the microscope: it will cause PERMANENT DAMAGE TO YOUR EYES!**
- Do not disassemble the camera on your own.
- Keep the camera away from moisture; do not use it in rainy weather.
- Keep the camera away from shock or excessive pressure.
- Do not overtighten the locking screws.
- Keep the camera away from hazardous environment, home and car heaters, incandescent lamps or open fire.
- When cleaning the lenses, first blow any dust or debris off the surface or wipe them off with a soft brush. Then wipe the lens with a soft cleaning cloth slightly damp with spirits or ether.

- Seek medical advice immediately if a small part or a battery is swallowed.

Battery safety instructions

- Always purchase the correct size and grade of battery most suitable for the intended use.
- Always replace the whole set of batteries at one time; taking care not to mix old and new ones, or batteries of different types.
- Clean the battery contacts and also those of the device prior to battery installation.
- Make sure the batteries are installed correctly with regard to polarity (+ and -).
- Remove batteries from equipment that is not to be used for an extended period of time.
- Remove used batteries promptly.
- Never attempt to recharge primary batteries as this may cause leakage, fire, or explosion.
- Never short-circuit batteries as this may lead to high temperatures, leakage, or explosion.
- Never heat batteries in order to revive them.
- Remember to switch off devices after use.
- Keep batteries out of the reach of children, to avoid risk of ingestion, suffocation, or poisoning.

Levenhuk Limited Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a lifetime warranty against defects in materials and workmanship. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for six months from date of retail purchase. Levenhuk will repair or replace such product or part thereof which, upon inspection by Levenhuk, is found to be defective in materials or workmanship. As a condition to the obligation of Levenhuk to repair or replace such product, the product must be returned to Levenhuk together with proof of purchase satisfactory to Levenhuk. This warranty does not cover consumable parts, such as batteries.

A Return Authorization (RA) Number must be obtained in advance of return. Contact the local Levenhuk branch to receive the RA number to be displayed on the outside of your shipping container. All returns must be accompanied by a written statement setting forth the name, address and telephone number of the owner, including a description of any claimed defects. Parts or products for which replacement is made will become the property of Levenhuk.

The customer will be responsible for all costs of transportation and insurance to and from Levenhuk or its authorized dealers and will be required to prepay such costs. Levenhuk will use reasonable efforts to repair or replace any product covered by this warranty within thirty days of receipt. If a repair or replacement will require more than thirty days, Levenhuk will notify the customer accordingly. Levenhuk reserves the right to replace any product that has been discontinued from its product line with a new product of comparable value and function.

This warranty does not apply to any defects or damages resulting from alteration, modification, neglect, misuse, usage of improper power sources, damage in transportation, abuse, or any cause other than normal use, or to malfunction

or deterioration due to normal wear. Levenhuk disclaims all warranties, express or implied, whether of merchantability or fitness for a particular use, except as expressly set forth herein. The sole obligation of Levenhuk under this limited warranty will be to repair or replace the covered product, in accordance with the terms set forth herein. Levenhuk disclaims liability for any loss of profits, loss of information, or for any general, special, direct, indirect or consequential damages which may result from breach of any warranty, or arising out of the use or inability to use any Levenhuk product. Any warranties which are implied and which cannot be disclaimed will be limited in duration to a term of six months for accessories from the date of retail purchase.

Some states/provinces do not allow the exclusion or limitation of incidental or consequential damages, so the above limitations and exclusions may not apply to you. This warranty gives you specific legal rights, and you may have other rights which vary from state to state or province to province.

Levenhuk reserves the right to modify or discontinue any product without prior notice.

NOTE: This warranty is valid to USA and Canadian customers who have purchased this product from an authorized Levenhuk dealer in the USA or Canada. Warranty outside the USA or Canada is valid only to customers who purchased from an authorized Levenhuk dealer in the specific country or international distributor. Please contact them for any warranty service.

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch:

Levenhuk Worldwide:

USA: www.levenhuk.com

Canada: www.levenhuk.ca

Czech Republic: www.levenhuk.cz

Netherlands: www.levenhuk.nl

Poland: www.levenhukoptics.pl

Russia: www.levenhuk.ru

Ukraine: www.levenhuk.ua

EU: www.levenhuk.eu

Purchase date _____ Signature _____ Stamp

Camera model	C35 NG	C130 NG	C310 NG	C510 NG	C800 NG	C1400 NG
Max. resolution (for still images)	640x80	1280x1008	2048x1536	2592x1944	3264x2448	4096x3288
Megapixels	0.3	1.3	3.0	5.0	8.0	14.0
Sensor	1/3" CMOS	1/2" CMOS	1/2" CMOS	1/2.2" CMOS	1/2.2" CMOS	1/2.3" CMOS
Pixel size, μm	5.6x5.6	3.6x3.6	3.2x3.2	2.2x2.2	1.67x1.67	1.4x1.4
Sensitivity, v / lux.sec @ 550 nm	1.9	1.0	1.0	0.53	0.3	0.724
Video frame rate, fps	<80	<50	<43	<60	<27	<27
Video recording	yes					
Dynamic range, dB	60	71	61	66.5	34	65.3
Mounting location	eyepiece tube (replaces the eyepiece)					
Image format	*.jpg, *.bmp, *.png, *.tif, *.gif, *.psd, *.ico etc.					
Spectral range, nm	380-650 (built-in infrared filter)					
Exposure	ERS (Electronic Rolling Shutter)					
White balance	auto/manual					
Exposure control	auto/manual					
Image control	image size, brightness, exposure control					
Interface	USB 2.0, 480Mb/s					
System requirements	Mac OS, Linux, Windows XP/Vista/7/8 (32-bit or 64-bit); USB port 2.0					
Software	USB 2.0 driver, Levenhuk ToupView software for image viewing and editing					
Casing	metal, painted black with a gold trim					
Power supply	via USB 2.0 cable					
Operating temperature range, °C	-10... 50					
°F	14... 122					

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Levenhuk® is a registered trademark of Levenhuk, Inc.
© 2006-2014 Levenhuk, Inc. All rights reserved.
Levenhuk, Inc. 1935 Brandon Court, Suite A-1